

## AntiMycoplasma Solution (Catalog#TBS8006)

### BACKGROUND

Mycoplasma contamination is one of the most common and major problems in cell culture for research laboratories and industries. A conservative estimate states that around 15-35% of all continuous cell cultures are contaminated with mycoplasma. Mycoplasma is one type of the smallest and simplest self-replicating prokaryotes, which can't be detected by visual inspection and gradually effect cell growth and alter DNA, RNA, and protein synthesis of host cells.

### DESCRIPTION

AntiMycoplasma Solution is designed to clean up mycoplasma contamination from cultured cells when it has been added into culture medium. The activity of solution is stable in culture medium and not affected by penicillin, streptomycin, and serum. This solution exhibits no toxicity to cultured cells.

### CONTENT

AntiMycoplasma Solution is supplied in 2x1ml tubes at a concentration of 25mg/ml.

### STERILITY

The solution is sterilized and filtered with 0.2 µm filter.

### APPLICATIONS

Treat mycoplasma contaminated cells.

### STORAGE CONDITIONS

AntiMycoplasma Solution is shipped at room temperature and should be stored at -20°C upon the arrival for long-term application. The solution is stable for 2 weeks at room temperature and at least one year at -20°C. Avoid repeated freeze-thaw cycles.

This product is for in vitro research use only and is not intended for use in humans or animals in therapeutic or diagnostic procedures.

### SUGGESTED PROTOCOL

To treat mycoplasma contaminated cells, the environment should be cleaned before treatment.

1. Decontaminate the external surface of the tube of solution by spraying with 70% ethanol.
2. Add AntiMycoplasma Solution into culture medium at a concentration of 25µg/ml (1:1000 dilution from stock concentration of 25mg/ml) or 12.5µg/ml (1:2000 dilution from stock concentration of 25mg/ml) under aseptic conditions.
3. Remove the old medium and replace with fresh medium containing AntiMycoplasma Solution as volume as needed.
4. The medium should be changed every other day. If the cells are stem cell or iPS cell, the medium should be changed every day.
5. The cells should be treated at concentration of 25µg/ml for 2 weeks or at concentration of 12.5µg/ml for 3 weeks.
6. The cells should be tested by using conventional PCR or cell based staining following the treatment.
7. To maintain the mycoplasma free culture, use Mycoplasma Prevention Solution (Cat# TBS8005) in the cell culture.

### REFERENCES

1. Drexler HG, Uphof CC (2002). Cytotechnology 39: 75-90
2. Aldecoa-Otalora E, Langdon WB, Cunningham P, Arno MJ (December 2009). "Unexpected presence of mycoplasma probes on human microarrays". BioTechniques 47 (6): 1013-5.

### RELATED PRODUCTS

Mycoplasma Prevention Solution(TBS8005)  
ESC/iPSC-qualified FBS (TBS8002)  
Fetal Bovine Serum (Cat#TBS8003)  
PBS (Cat#TBS5003, -01, -02or TBS5027)

#### Contact

Phone: 650-917-9269

Email: [info@tribioscience.com](mailto:info@tribioscience.com)

Fax: 650-618-5498

Web: [www.tribioscience.com](http://www.tribioscience.com)