

## Probe qPCR Detecting GMO Marker 98140 from Maize

### DESCRIPTION

Many countries in the world regulate the cultivation and trade with genetically modified organisms (GMO). The enforcement of such regulations depends on the ability to detect and quantify the presence of GMOs in food, feed, and seed products. Real-time qPCR method is sensitive and robust, and is regarded as the gold standard for GMO analysis.

The Maize GMO 98140 qPCR Kit is designed for identifying GMO marker 98140 presence or absence in food, feed and seed products using real-time quantitative polymerase chain reaction (qPCR), primers, and labeled probe. The Kit includes maize GMO 98140 Positive and Negative controls, PCR internal controls, qPCR Super Mix (M1), maize GMO 98140 Prime-Probe Mix (M2), in which the probe is labeled with Fam, and Hex is labeled for PCR internal control. These aids in the straightforward interpretation of the results.

### PRINCIPLE

GMO marker identification is based on qPCR amplification of a specific region of the relevant target genome. The amplified product is detected using target-specific fluorescent probes that bind to the amplified product. As the PCR product accumulates, there is an increased fluorescence signal from the bound probes. Monitoring the fluorescence intensities during the PCR run allows the detection of the accumulating PCR product in real time.

### KEY FEATURES

- ❖ Highly sensitivity and specificity for identification of GMO 98140.
- ❖ High efficiency: the optimal systemic conditions for PCR amplifications.
- ❖ Streamlined protocol: Just add DNA Template, and water.
- ❖ No cross reactivity with others.

### APPLICATIONS

Detect GMO marker in food, and animal feed.

### KIT CONTENTS

Name	100RXN
qPCP Super Mix	0.8 mL
Primer-probe Mix	0.6 mL
Positive Control DNA	60 µL
Negative Control DNA	60 µL

GMO98140 gene is labeled with FAM;  
PCR internal control is labeled with Hex.

### STORAGE CONDITION

The kit is shipped on ice and stored at -20°C for long-term storage. Shelf life of 12 months after receipt.

### PCR PROTOCOL

1. **DNA extraction:** The methods for DNA extraction can be used for any suitable preparation of DNA purification from food samples. We recommend that the Fast genomic DNA extraction method is used for this purpose (catalog: TBS6008).
2. **Set up a PCR reaction for each sample in 20 µL**

Reaction Component	Volume (µL)
qPCR Super Mix	7.0
Primer-probe Mix	5.0
Nuclease-free Water	3.0
DNA sample	5.0
<b>Final Volume</b>	<b>20 µL</b>

Internal control should be included as below: Positive Control or Negative control (5 µL /reaction).

### 3. Suggested PCR conditions

Step	Amplification	PCR	
	HOLD	CYCLE (40 cycles)	
		Denature	Anneal/ Extend
<b>Temperature</b>	95 °C	95 °C	60 °C
<b>Time</b>	1 min	15 sec	60 sec

### DATA ANALYSIS

Positive Reaction: Sample Ct < or = 37, and Positive, Negative and Blank controls are normal.

Negative Reaction: Sample Ct ≥ 38, and Positive, Negative and Blank controls are normal.

PCR internal control is positive in all samples, positive and negative controls. The positive response indicates a normal PCR amplification. Otherwise, the PCR reaction may be inhibited.

Repeat Reaction: If one of the control reactions is not normal, PCR reaction is failed, and should be repeated.

### RELATIVE PRODUCTS

- TBS6008: Fast Genomic DNA Extraction
- TBS6025: Microbial DNA Magnetic Extraction
- TBS43002: Maize GMO CBH-351 qPCR Detection
- TBS43003: Maize GMO DBT418 qPCR Detection
- TBS43004: Maize GMO LY038 qPCR Detection
- TBS43005: Maize GMO Mon863 qPCR Detection
- TBS43006: Maize GMO Mon87411 qPCR Detection
- TBS43007: DAS-59132 qPCR Detection
- TBS42025: 4-In-1 Aspergillus Species qPCR
- TBS42026: O157H7 E. Coli qPCR
- TBS42028: Salmonella qPCR
- TBS42029: Salmonella-STEC qPCR
- TBS42031: Listeria Monocytogen qPCR
- TBS42032: Listeria Species qPCR
- TBS42033: Bacillus Cereus qPCR

**For research use only.**